



MagReSyn[®] Streptavidin MS: enabling more sensitive affinity workflows

Aaron Fehr, Emanuela Milani, Bernd Wollscheid
Institute of Translational Medicine, Dept of Health Sciences
and Technology, ETH Zurich

Isak Gerber, Stoyan Stoychev, Andrea Ellero, Justin Jordaen
ReSyn Biosciences (Pty) Ltd

HIGHLIGHTS

- Protease-resistant MagReSyn[®] Streptavidin MS engineered for applications requiring on-bead digestion
- Multiple-fold reduction in streptavidin contaminating peptides during on-bead trypsinization
- Fewer contaminating peptides compact dynamic range thereby improving detection of target peptides
- Reduced non-specific binding of background proteins and peptides compared to standard product
- Retention of >80% of the capacity of standard MagReSyn[®] Streptavidin
- Workflow is fully automated on the Agilent Bravo liquid handling system


INTRODUCTION

To understand cellular function, it is critical to characterize interactions among biomolecules. One common strategy for proximity-based protein interaction analysis (APEX, BioID, LUX-MS etc) is based on the enrichment of biotinylated peptides/proteins using a streptavidin matrix, such as MagReSyn[®] Streptavidin. The very high affinity between biotin-tagged targets and streptavidin allows for the enrichment and subsequent detection of interaction partners while stringent washing conditions reduce non-specific binding. This strategy works well if the bound proteins/biomolecules can be eluted using harsh conditions, such as boiling the MagReSyn[®] Streptavidin beads in Laemmli buffer.

However, for mass-spectrometry-based (MS) identification, an on-bead digestion of bead-bound biotinylated proteins is typically preferred/employed. In such a digestion process, affinity-enriched proteins are digested, but also parts of the streptavidin matrix. In the subsequent LC-MS analysis, streptavidin-derived peptides mask the detection of low-abundance peptides of interest.

To circumvent these concerns, the new MagReSyn[®] Streptavidin MS is specifically designed to cater to applications requiring on-bead enzymatic digestion of proteins. This is achieved through the chemical modification of lysine and arginine residues of the immobilized Streptavidin as described by Rafiee *et al.* in 2020. (DOI: [10.15252/msb.20199370](https://doi.org/10.15252/msb.20199370)). The derivatization was optimized to retain a high affinity for biotin, retaining over 80% of the capacity of the standard MagReSyn[®] Streptavidin product, with multiple-fold decreases in the release of streptavidin peptides resulting from on-bead digestion.

METHODS

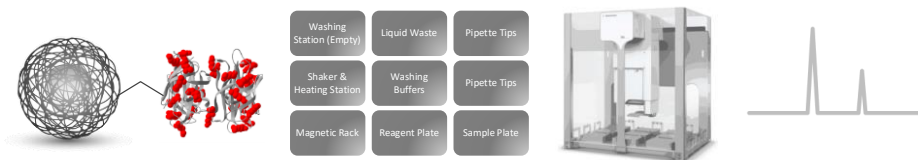

Evaluating binding capacity and reduction in streptavidin peptides from on-bead digest



- The binding capacity of MagReSyn® Streptavidin MS was determined by incubating beads with pure biotinylated rabbit IgG and biotinylated oligo (15mer) to estimate capacity before and after chemical treatment of lysine and arginine residues.
- On-bead digestion was performed by incubating 20 µl of MagReSyn® Streptavidin or MagReSyn® Streptavidin MS, pre-equilibrated in 25 mM Ammonium Bicarbonate pH 8, with Trypsin (1:20) for 16 hours at 37°C.
- Digests were loaded on 75 µm x 2 cm Acclaim™ PepMap™ trap column using 2% ACN, 0.2% FA and separated via 15 min linear gradient on Waters CSH 75 µm x 25 cm column coupled to Sciex 6600 TToF operated in SWATH mode (48VW).
- Targeted data extraction for five streptavidin-derived peptides was performed in Skyline with sum of transition peak area utilized as a measure of peptide abundance.

Binding capacity and reduction in streptavidin-derived peptides in a spiked cellular lysate


- The workflow was automated on an Agilent Bravo liquid handling station.
- Biotinylated target molecules were captured by incubating 15 µl MagReSyn® Streptavidin, or 16.5 µl MagReSyn® Streptavidin MS (corrected for capacity) with 400 µg of 1% NHS-biotinylated yeast protein spiked into HEK lysate for 20 minutes (quadruplicate samples).
- Non-specifically bound proteins were removed by 3 consecutive washes consisting of:
 - Wash 1: 1% SDS in PBS
 - Wash 2: 8M Urea in 50 mM Ammonium Bicarbonate (pH 8).
 - Wash 3: 5 M NaCl
 - Wash 4: 100 mM NaHCO₃
 - Wash 5: 80% Isopropanol
 - Wash 6: 100 mM Ammonium Bicarbonate (pH 8).
- Target release was achieved via on-bead digestion performed overnight in 100 µl of 0.1% RapiGest™ in 100 mM Ammonium Bicarbonate digest buffer using 0.8 µg of Trypsin.
- Peptides were desalted using C18 cartridges, dried and resuspended in 3% ACN, 0.1% FA. Equal amounts were separated at 250 nl/min over a 70 min gradient of 5 to 41% ACN, 0.1% formic acid (FA), on a 75 µm x 15 cm, in-house packed C18-column connected to a nano-flow HPLC (ThermoFisher Scientific). Mass spectra were acquired on an Orbitrap Q Exactive™+ operated in DIA mode using 18 m/z windows and staggered collision energy (nce=25,27,30).
- Raw data files were processed using DIA-NN with an in-silico-generated spectral library. All reported data was filtered at 1% FDR and normalized intensities are shown.

RESULTS

Binding Capacity: MagReSyn® Streptavidin MS showed a marginal decrease in binding capacity of <15% in comparison to the standard Streptavidin (below left), but with a concomitant >95% reduction in abundance of streptavidin-derived peptides after on-bead digestion (below right).

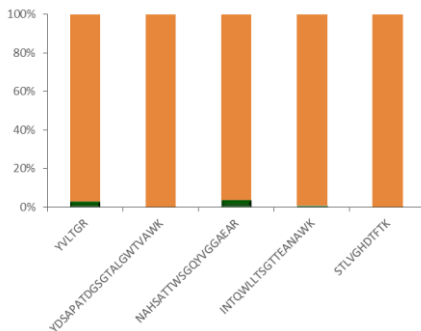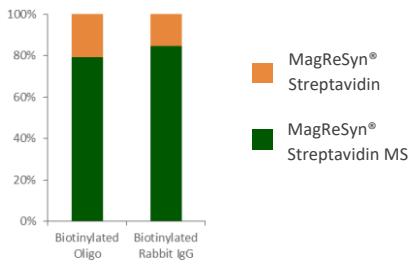




Summed Precursor Quantities

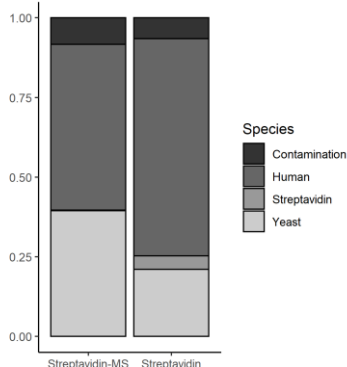

Isolation of biotinylated yeast proteins: Fewer streptavidin precursors, and with significantly lower intensity, were detected after the on-bead digestion with MagReSyn® Streptavidin MS after isolating biotinylated yeast proteins from a complex HEK cell lysate. This resulted in a higher number of yeast-derived target precursors, while the number of contaminating human background precursors decreased.

Samples enriched using MagReSyn® Streptavidin MS resulted in a similar number of yeast proteins being identified, but with more peptides isolated per protein, while the number of human-derived, contaminant peptides decreased significantly.

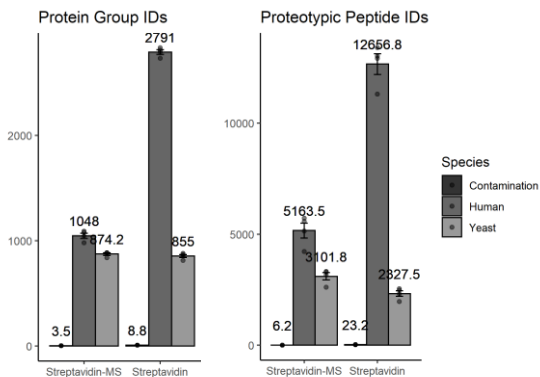

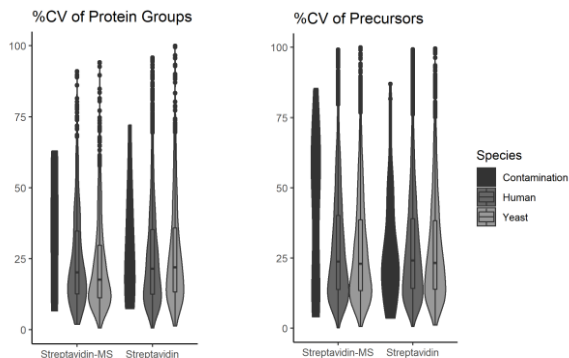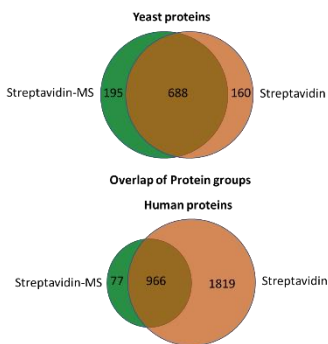

Over 80% of the target proteins detected in yeast overlapped between the standard Streptavidin and Streptavidin MS products (above left), but the technical reproducibility increased with median CV of 18% (Streptavidin MS) versus 22 % (standard Streptavidin – violin plots above).

CONCLUSIONS

- ✓ We describe a new digestion-resistant MagReSyn® Streptavidin MS product
- ✓ The binding capacity for biotinylated IgG shows only a marginal decrease, but with the vast reduction in the abundance of streptavidin-derived peptides after on-bead digestion
- ✓ When evaluated for the capture of NHS-biotinylated yeast proteins in a complex lysate, the result was an increased number of identifications and increased confidence in quantitation.
- ✓ This in turn resulted in an overall increase in the reproducibility of the workflow.
- ✓ The chemical modification of lysine and arginine residues in streptavidin, had a further positive effect by improving the selectivity over non-treated, which could possibly be attributed to reduction in ionic interactions with the lysate.

REFERENCES

R Studio – Available online: <https://posit.co/download/rstudio-desktop/>
 Rafiee et al., 2020. Mol. Systems Biol. DOI: 10.15252/msb.20199370

MagReSyn is a registered trademark of ReSyn Biosciences (Pty) Ltd, Acclaim, Pepmap & Q Exactive are trademarks of Thermo Fisher Scientific, Inc. and its subsidiaries. RapiGest is a trademark of Waters Corporation,

Agilent Bravo scripts available on request.